

# NOCOBOX



Manufactured in France by Oxy'Pharm according to ISO 13485.

The **Nocobox** are culture dishes with nutrients to highlight germs on surfaces. They are used to verify the result of disinfection from an airborne surface disinfection system. They contain a nutrient that promotes the growth of bacteria and therefore highlights their presence.


## STORAGE

Store the boxes in a cool (below 30°C) and dry place. Store out of sunlight. Length of validity in the original closed packaging: 4 months.

Before



After

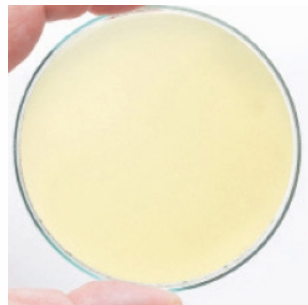

## IMPORTANT

Elimination: collection and incineration by specialised companies.

## INSTRUCTIONS FOR USE

### 1<sup>st</sup> sampling before disinfection

- 1 Open the box carefully (do not touch the test area).
- 2 Apply 10 seconds of light pressure on the surface to be tested.
- 3 Close the box.
- 4 Leave it at development temperature (20°C - 25°C) for 48 hours.
- 5 Read the results.

### 2<sup>nd</sup> sampling after disinfection

- 6 Repeat steps 1 to 5 after disinfection using new boxes.
- 7 Compare and observe the effectiveness of the treatment.

## REFERENCES

### PACKAGING

### REFERENCE

Boxes of 10

5000.020